K-SYSTEMS

USER MANUAL

РУКОВОДСТВО ПОЛЬЗОВАТЕЛЯ (РУССКИЙ, RUSSIAN)
BENUTZERHANDBUCH (DEUTSCH/GERMAN)
MANUAL DEL USUARIO (ESPAÑOL/SPANISH)
MODE D'EMPLOI (FRANÇAIS/FRENCH)
MANUALE UTENTE (ITALIANO/ITALIAN)
用户手册 (简体中文 / CHINESE)

Anti Vibration Table - AV1 and AV5 (English)

Антивибрационный стол AV1 и AV5 (РУССКИЙ, RUSSIAN)
Antivibrationstisch – AV1 und AV5 (DEUTSCH/GERMAN)
Mesa antivibración - AV1 y AV5 (ESPAÑOL/SPANISH)
Table antivibration - AV1 et AV5 (FRANÇAIS/FRENCH)
Tavolo antivibrazione: AV1 e AV5 (ITALIANO/ITALIAN)
防振动工作台 - AV1 和 AV5 (简体中文 / CHINESE)

AV1 AND AV5 USER MANUAL

AV1 AND AV5 USER MANUAL (English).....	3
РУКОВОДСТВО ПОЛЬЗОВАТЕЛЯ AV1 И AV5 (РУССКИЙ, RUSSIAN).....	14
BENUTZERHANDBÜCHER FÜR AV1 UND AV5 (DEUTSCH/GERMAN)	25
MANUAL DEL USUARIO DE AV1 Y AV5 (ESPAÑOL/SPANISH).....	36
MODE D'EMPLOI DES TABLES AV1 ET AV5 (FRANÇAIS/FRENCH)	47
MANUALE UTENTE AV1 E AV5 (ITALIANO/ITALIAN).....	58
AV1 和 AV5 用户手册 (简体中文 / CHINESE)	69

CONTENT

1. GENERAL INFORMATION AND SERVICE	4
1.1 General Description	4
1.2 Copyright	4
1.3 Customer Service Contact Information	4
2. SETUP	5
2.1 Unpacking and Inspection	5
2.2 Assembling the AV1/AV5	5
2.3 Weight Adjustment	7
2.3.1 Removing the Middle Spring	8
2.3.2 Adding the Weight Plate	9
3. TROUBLESHOOTING	10
4. MAINTENANCE	11
5. TECHNICAL DATA AV1/AV5	11
6. WARRANTY INFORMATION AND LIMITS ON LIABILITY	12

1. GENERAL INFORMATION AND SERVICE

1.1 General Description

The K-SYSTEMS Anti Vibration tables are designed to meet the requirements of ICSI Procedures. With an isolated microscope platform and a complex anti-vibration system, the table is preventing vibrations from reaching the microscope platform and dampening unavoidable vibrations in the shortest possible time.

The inverted microscope is placed on an individually suspended floating platform. The platform must be levelled and isolated – meaning not touching – the rest of the table plate. The vibrations are reduced by friction between the platform and the springs holding the platform.

1.2 Copyright

This manual contains information that is subject to copyright. All rights reserved. This manual should not be photocopied, otherwise copied or distributed, completely or in part, without the approval of CooperSurgical, Inc.

Users of K-SYSTEMS products should not hesitate to contact us if there are any unclear points or ambiguities in this manual.

1.3 Customer Service Contact Information

European Service Representatives

ORIGIO a/s
Knardrupvej 2
2760 Måløv
Denmark

Customer Service:
+45 46 79 02 02
Fax: +45 46 79 03 02

Email:
New Purchases:
customerservice@origio.com

Service Related Emails:

service@origio.com
www.origio.com

U.S. Service Representatives

CooperSurgical, Inc.
95 Corporate Drive Trumbull, CT
06611 USA

Customer Service:
Phone: (800) 243-2974
Fax: (800) 262-0105

International
Phone: +1 (203) 601-9818
Fax: +1 (203) 601-4747

Service Related Emails:

service@origio.us.com
www.coopersurgical.com

2. SETUP

2.1 Unpacking and Inspection

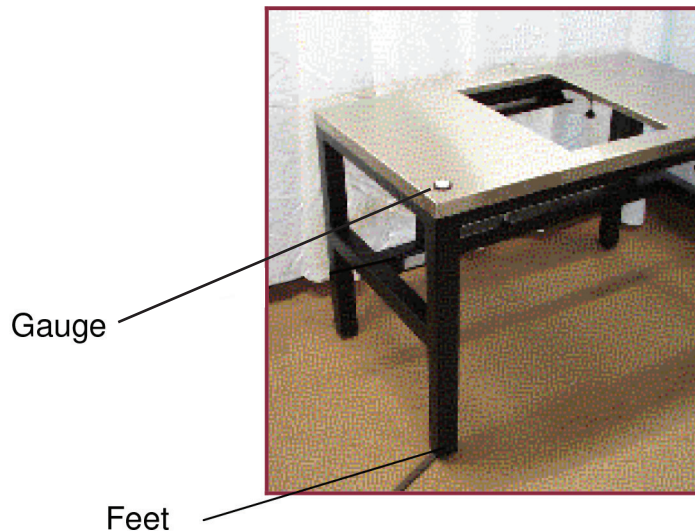
- Unpack the equipment.
- Check the packing and picking list to insure all accessories are there.
- Examine the packing material carefully for separately packed items.
- Check the table for external surface damage.
- Please be aware that the springs must not be touched at any time. Please use gloves when you assemble the unit.

2.2 Assembling the AV1/AV5

For assembling the AV1 /AV5 please avoid touching the springs.

Place the anti vibration table in its required position on the floor. Preferably on a hard floor.

Level the table plate horizontally by adjusting the feet on the table legs. To ensure the best result, place a gauge/bulls eye on the table plate and make sure it is leveled.

NOTE: The table can be modified to optimally work with equipment of different weight. It is important that you know the exact weight of the equipment before proceeding. **See chapter 3.3.**

Put on the gloves. Remove the foam pads from the springs – without touching the springs with your bare hands.

Place the floating plate upon the springs, leave equal distance from all sides to the hole of the table plate for centering the floating plate. It is very important that there is no contact between the floating plate and the table!

Place the inverted microscope on the floating plate.

The position and height of the floating plate are adjusted by:

Loosen the four securing bolts

Turn the adjustment screws to the necessary position.

Tighten the four securing bolts again.

2.3 Weight Adjustment

The anti-vibration table AV1 and AV5 can be used to hold equipment with a weight from 15 to 72 kg. However, the table needs to be installed correctly to have the optimal anti vibration ability.

NOTE: Before proceeding it is important that you know the exact weight (in Kg) of the equipment, which you will place on the anti-vibration platform.

When AV1 or AV5 leaves the factory it is equipped with 3 springs and an extra weight plate. See right column in the table below. The yellow areas indicate the optimal working conditions for the 3 different installation methods.

Example: The total weight of equipment placed on the anti-vibration platform with 3 springs should be in the range from 45- 72 Kg.

The Anti-vibration table can be modified in 2 steps to gradually work optimal with less heavy equipment on the platform.

Step 1: To work with equipment from 30-45 kg one spring is removed so only 2 springs are used.

NOTE: Only remove the middle spring, when needed.

Step 2: To work optimal with equipment from 15-30 kg 2 springs are installed and an extra weight plate must be used.

		2 Springs and Weight Plate		2 Springs without Weight Plate		3 Springs without Weight Plate (From Factory)	
Equipment weight	Use:	Deflection	Frequency	Deflection	Frequency	Deflection	Frequency
Kg		mm	Hz	mm	Hz	mm	Hz
15-20	2 Springs and Weight Plate	15.59	3.99				
20-25		18.13	3.70				
25-30		19.94	3.53	15.59	3.99	10.40	4.89
30-35	2 Springs without Weight Plate	21.76	3.38	18.13	3.70	12.09	4.54
35-40		23.57	3.25	19.94	3.53	13.30	4.33
40-45				21.76	3.38	14.51	4.14
45-50	3 Springs without Weight Plate			23.57	3.25	15.71	3.98
50-55				25.38	3.13	16.92	3.83
55-60						18.13	3.70
60-65						19.34	3.59
65-70						20.55	3.48
70-72						21.76	3.38

Instructions about how to remove the middle spring and how to install the extra weight plate can be found in the following pages:

2.3.1 Removing the Middle Spring

If the weight of the equipment that the AV table is going to hold is between 27 and 47 kg the middle spring needs to be removed. This is done in the following manner:

1. Unscrew the table plate; it is attached with 4 screws on the bottom side (only on AV1).
2. Remove the table plate (only on AV1).
3. Please put on gloves when touching the springs
4. In each side of the table there are metal rods screwed into the frame, unscrew these two metal rods.

5. Push the springs to one end and lift out the other metal rod

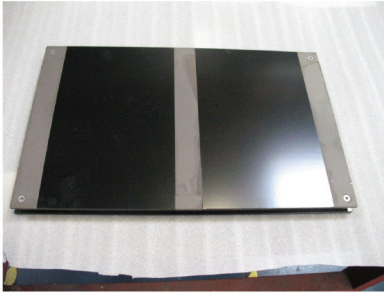
6. Remove the screw in the middle spring on the opposite side of the end that is lifted.
7. Pull out the middle spring.
8. Push the two remaining springs to the bottom side.
9. Move the other metal rod in place and screw both metal rods in place again.
10. Place the table plate back on and attach it underneath with the 4 screws.

2.3.2 Adding the Weight Plate

AV1 Table


- 1) Place the floater on a flat surface


- 2) Remove the screws from the bottom plate, remove the plate.

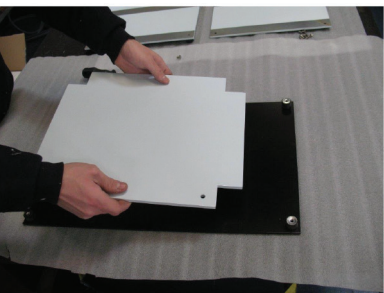- 3) Add the extra weight

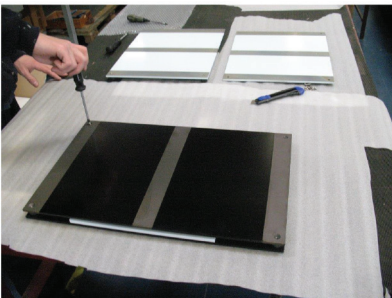
- 4) Mount and screw the bottom plate on again


AV5 Table

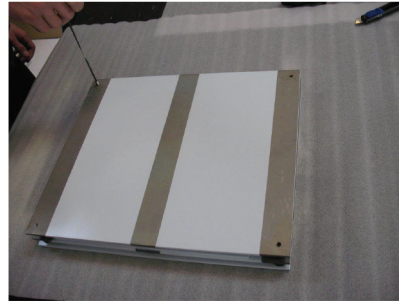
- 1) Place the floater on a flat surface and remove the screws

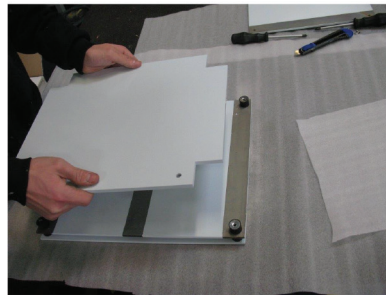
- 2) Remove the bottom plate and add the extra weight

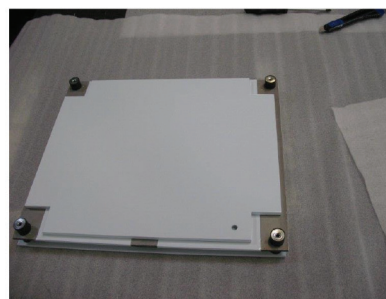- 3) The extra weight in place

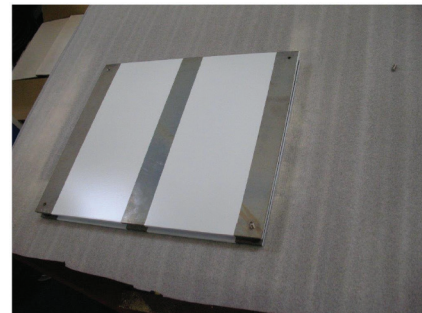- 4) Mount and screw the bottom plate and reinsert screws

3. TROUBLESHOOTING

VIBRATIONS CAUSED BY	ACTION TO BE TAKEN
The floater is touching the rest of the table plate and vibrations are being transferred directly to the microscope.	Center the floater within the hole. Air should be all around the floater.
Loose tightening bolts.	Tighten the four tightening bolts.
The spring system is not painted – but need to be spotless clean and so should the floater where it rests on the springs. If there is dirt on any of these areas, friction is lost and vibrations are transferred to the microscope.	Clean the areas with 70% alcohol.
Tight cables from the micromanipulators will cause vibrations to be transmitted directly from the micromanipulator to the microscope.	Loosen the cables so they hang loose.
The floater is placed too high, resulting in the springs touching the underside of the table plate.	Lower the floater so it is in level with the rest of the table plate.
Microscope not placed evenly on the floater.	The weight of the microscope needs to be distributed evenly on the floater. Since most microscopes are heaviest in the back part, the microscope should be moved more to the front of the table plate.

4. MAINTENANCE

All surfaces should be cleaned with a 70% alcohol solution on a clean cloth or lintfree paper towel.

1. Clean all components by wiping gently with gauze. To remove fingerprints or oil smudges, wipe with gauze slightly moistened with a mixture of ether (70%) and alcohol (30%).

Since solvents such as alcohol are highly flammable, it must be handled carefully. Be sure to keep these chemicals away from open flames or potential sources of ignition – for example, electrical equipment that is being switched on or off. Also remember to always use these chemicals only in a well ventilated room.

2. If smudges are difficult to clean, wipe them with a soft cloth slightly moistened with a diluted neutral detergent.

5. TECHNICAL DATA AV1/AV5

Material.....	Stainless steel
Weight AV1/AV5.....	100Kg
Stainless steel plate AV1/AV5.....	1200x800 mm
Height of table plate	780 mm
Undamped frequency	< 4 Hz
Maximum deflection.....	22 mm
Mass (to be dampened)	15-72 Kg

6. WARRANTY INFORMATION AND LIMITS ON LIABILITY

CooperSurgical warrants that the workstation will be free from defects in materials and workmanship for one year from the date of initial purchase.

If CooperSurgical determines that a workstation fails to conform to that warranty during the one-year warranty period, CooperSurgical will, as the sole remedy for that failure to conform, repair or replace that workstation, at CooperSurgical's discretion, free of charge.

To return a workstation to CooperSurgical, a customer must comply with CooperSurgical's Returned Goods Policy described in this manual. A customer will not have any remedy if a workstation does not conform to the warranty for that workstation unless that workstation is returned to CooperSurgical in accordance with that Returned Goods Policy.

CooperSurgical will ship returned products that it repaired or replaced under warranty to the customer who returned those products, at CooperSurgical's expense F.O.B. the customer's facility. Under all other circumstances, CooperSurgical will ship returned products to the customer who returned those products at the customer's expense F.O.B. CooperSurgical's facility.

CooperSurgical's warranties do not cover damage caused by misuse, improper care, improper use of chemicals or cleaning methods, loss, theft, servicing by non-authorized personnel or negligent or intentional conduct on the part of the owner or user of the workstation, nor do they cover normal wear and tear or general maintenance. Any modifications or changes to a workstation will void that workstation's warranty. CooperSurgical's warranties do not apply to any single- or limited-use, disposable or consumable components or items.

CooperSurgical is not responsible for, and the owner and operator of the workstation shall defend, indemnify and hold harmless CooperSurgical from and against, all claims, damages, and other losses resulting from the improper servicing, maintenance, repair use or operation of the workstation or the owner or operator's negligence or willful misconduct.

THE ABOVE WARRANTIES ARE IN LIEU OF, AND COOPERSURGICAL HEREBY DISCLAIMS, ALL OTHER WARRANTIES, EXPRESS OR IMPLIED, WRITTEN OR ORAL, WITH RESPECT TO COOPERSURGICAL'S PRODUCTS, INCLUDING THE WARRANTIES OF MERCHANTABILITY AND FITNESS FOR A PARTICULAR PURPOSE. NO TERMS, CONDITIONS, UNDERSTANDINGS OR AGREEMENTS THAT PURPORT TO MODIFY THE ABOVE WARRANTIES OR THAT MAKE ANY ADDITIONAL WARRANTIES FOR ANY COOPERSURGICAL PRODUCT SHALL HAVE ANY LEGAL EFFECT UNLESS MADE IN WRITING AND SIGNED BY AN AUTHORIZED COOPERSURGICAL CORPORATE OFFICER.

COOPERSURGICAL SHALL NOT UNDER ANY CIRCUMSTANCES BE LIABLE FOR LOST PROFITS, DAMAGES FROM LOSS OF USE OR LOST DATA, OR INDIRECT, SPECIAL, INCIDENTAL OR CONSEQUENTIAL DAMAGES UNDER ITS WARRANTIES OR OTHERWISE FOR ANY CLAIM RELATED TO COOPERSURGICAL'S PRODUCTS, EVEN IF COOPERSURGICAL HAS BEEN ADVISED, KNEW OR SHOULD HAVE KNOWN OF THE POSSIBILITY OF SUCH DAMAGES. COOPERSURGICAL'S LIABILITY WITH RESPECT TO A PRODUCT COVERED BY A WARRANTY OR OTHERWISE SHALL BE LIMITED IN ALL CIRCUMSTANCES TO THE PURCHASE PRICE OF THAT PRODUCT.

Returned Goods Policy

1. Goods will be accepted for return for the following reasons:

- If shipment was made without the customer's authorization or order
- If incorrect items were shipped
- If defective items were shipped
- If defective goods are covered by the standard warranty

WARRANTY INFORMATION AND LIMITS ON LIABILITY CONTINUED

- 2. To return goods, you must contact a Customer Service Representative by telephone (+45 46 79 02 02) for a Returned Merchandise Authorization (RMA) number. Items will not be accepted without an RMA number.**

Please have the following information:

- Reason you wish to return the goods
- Quantity, description, part number, serial number of the goods
- Date of receipt of order
- Customer's purchase order and the CooperSurgical invoice number

All used goods must be cleaned and sterilized prior to shipment.

- 3. Shipment must be sent prepaid by the customer. Freight collect shipments will not be accepted, and goods will be returned to sender.**

- 4. If Customer intends to return equipment ordered in error, the following restocking charges and terms will apply:**

- 25 percent within 60 days from date of shipment
- Goods must be returned unused, in the original carton, and in marketable condition
- Refurbishing and replacement charges will be added to the restocking charges for damaged or missing items
- No return after 60 days
- No refund on sterile, single-use disposable products

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**CooperSurgical European Distribution Center B.V.